

To boldly flow with Qdot® nanocrystals

Now available: Qdot® nanocrystal–conjugated primary antibodies

Researchers today are trying to maximize information from their flow cytometry experiments by looking at more parameters in one sample. Qdot® nanocrystals provide a powerful way to increase your fluor selection from your available excitation sources. Qdot® nanocrystal conjugates will allow you to add 1–6 colors to your data acquisition from a violet laser and, in addition, provide the advantages of brightness and photostability.

The fluorescence properties of Qdot® nanocrystals are different from those of typical dye molecules. Typical fluorescent dyes have excitation and emission spectra with relatively small Stokes shifts, which means that the optimal excitation wavelength is close to the emission peak. Qdot® nanocrystals have broad absorbance spectra and are optimally excited by a UV or violet (405–407 nm) laser, although usable excitation can also be obtained with other sources such as the 488 nm laser. Their emission peaks are narrow and symmetrical, and do not change with excitation (Figure 1). Because of their spectral properties, Qdot® nanocrystals are brighter than most common fluor.

Qdot® nanocrystal conjugates may be used in the same way as conventional conjugates. Because staining conditions may vary, reagents should be titrated with samples to obtain optimal staining concentrations. Figure 2 shows typical profiles of human peripheral blood leukocytes (PBLs) stained with Qdot® nanocrystal conjugates specific for CD3, CD4, CD8, and CD14 antigens. The two-color combination was analyzed with <4% compensation between channels. In addition, Qdot® nanocrystals are compatible with common lysing, fixation, and permeabilization reagents, such as Cal-Lyse™ and FIX & PERM®.

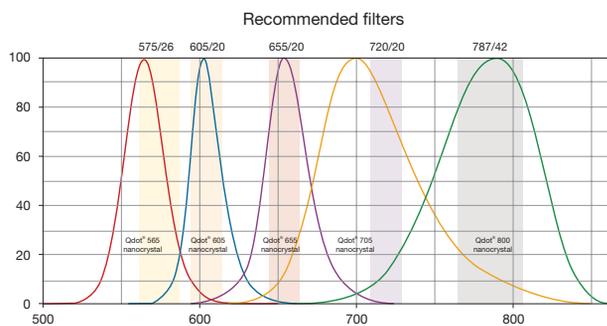


Figure 1—Recommended filter configuration and emission profiles for selected Qdot® nanocrystals. Filter diagrams and emission curves were viewed with the Fluorescence Spectra Viewer (probes.invitrogen.com/resources/spectraviewer/).

Primary antibody conjugates now available

As the exclusive provider of Qdot® nanocrystal technology for life science research, Invitrogen offers a full range of tools, from new primary antibody conjugates to secondary detection reagents, to maximize the use of your flow cytometer by combining Qdot® nanocrystal technology with existing organic fluorophores. For more information on using this new technology in flow cytometry, including detailed protocols and resources, visit www.invitrogen.com/qdotforflow.



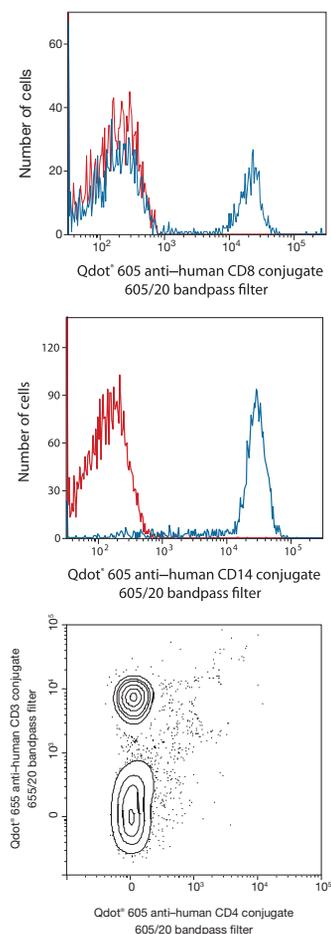


Figure 2—Staining profiles for Qdot® nanocrystal–conjugated antibodies. Human peripheral blood leukocytes were stained with the specified Qdot® conjugates. Samples were analyzed using a BD™ LSR II flow cytometer with 405 nm excitation and the specified emission filters.

Automated Affinity Purification

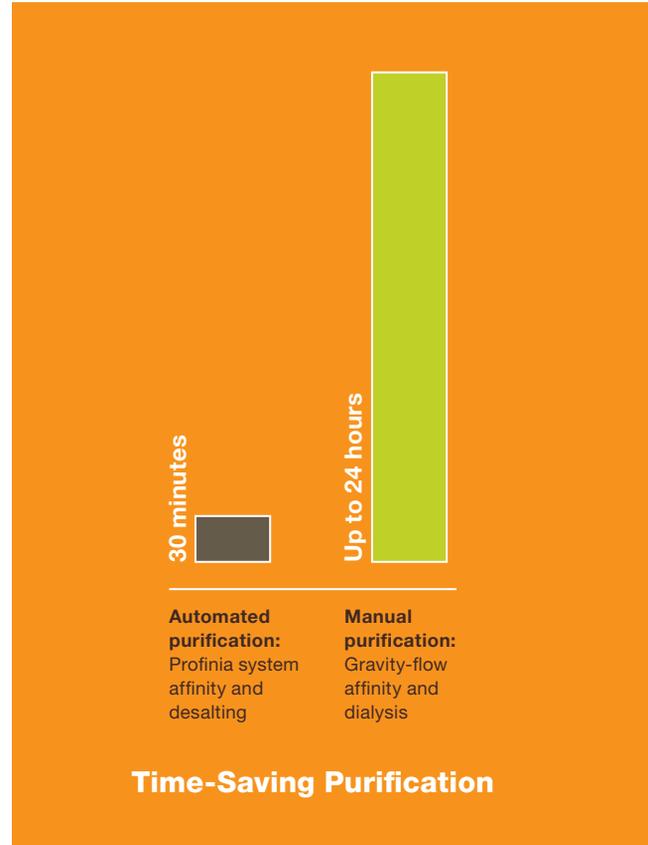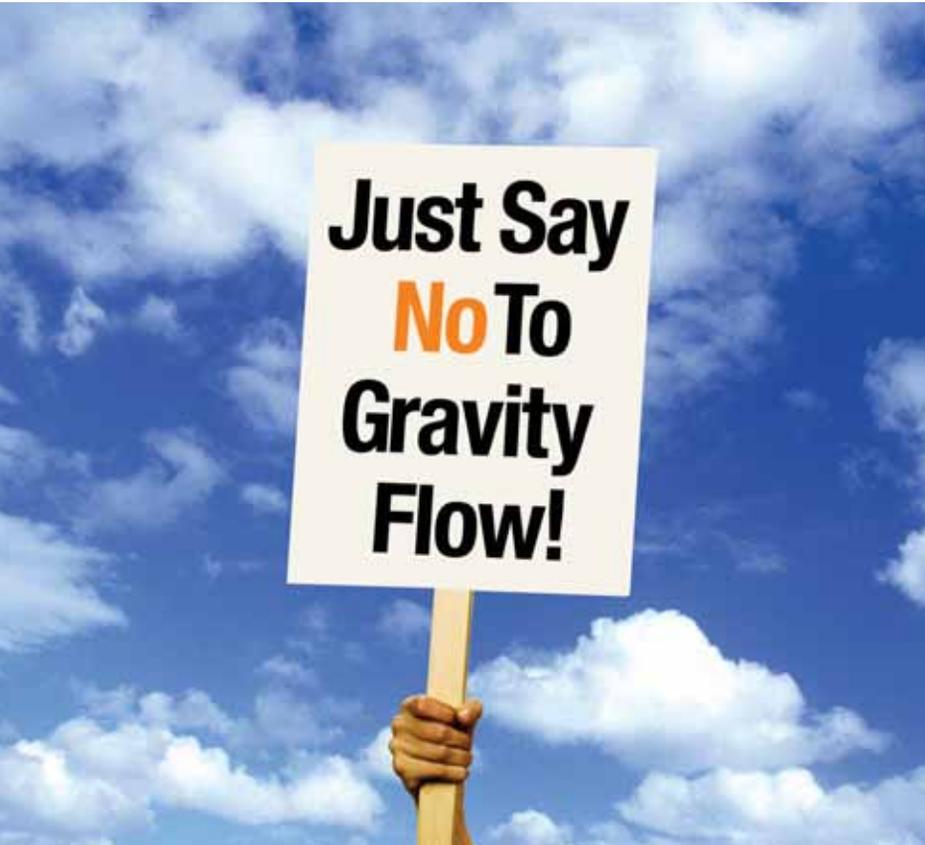

Say Yes to Automated Purification!

Free up hours of your day with the automated Profinia™ protein purification system.

A convenient and automated alternative to existing methods of purification, the Profinia protein purification system brings unprecedented speed and simplicity to the purification of affinity-tagged proteins.

- Preprogrammed methods for IMAC (native and denaturing), GST, and Protein A/G affinity purification, with integrated desalting applications
- Large, informative touch screen interface allows easy navigation through protocol steps
- Optimized kits and reagents match the methods and the instrument for greater reproducibility and reduced preparation time
- Automatically calculated run data include protein yield and concentration

For more information and a free IMAC, GST, or Protein A cartridge, visit us on the Web at www.bio-rad.com/ad/profinia/



Bio-Rad Laboratories, Inc.
Recipient of the 2007 Product Line Strategy Award,
U.S. Liquid Chromatography Market

Purification of fusion proteins may require a license from third parties.

To find your local sales office, visit www.bio-rad.com/contact/
In the U.S., call toll free at 1-800-4BIORAD (1-800-424-6723)

Visit us on the Web at discover.bio-rad.com



Now **more** than ever.



MoFlo™ XDP Cell Sorter



CyAn™ ADP Analyzer



FC 500 MPL Analyzer



EPICS™ XL-MCL™ Analyzer



Quanta™ SC MPL Analyzer



Vi-CELL® Cell Viability Analyzer



Z™ Series Coulter Counter



Think Beckman Coulter for flow cytometry.

No one can meet your diverse cell analysis needs like Beckman Coulter. For flexibility and reliability, depend on our expanding family of flow cytometry solutions that range from affordable analyzers to a cell sorter that sets the standard. You can always count on us—whether you're looking for instrumentation, reagents for your research and clinical needs, or the security of Beckman Coulter's global network of service and support. Discover all you need all in one place at www.beckmancoulter.com/flowcytometry



Genomics Proteomics **Cell Analysis** Centrifugation Lab Tools Particle Characterization Bioseparation Lab Automation

Simplify • Automate • Innovate

iSBTc

23rd Annual Meeting & Associated Programs

October 29 - November 2, 2008
San Diego, CA

www.isbtc.org

INTERNATIONAL SOCIETY FOR BIOLOGICAL THERAPY OF CANCER

**Abstract
Deadline:**

July 21, 2008

Program/Abstract Topics

- Adoptive Transfer
- Cancer and Inflammation
- Cancer Stem Cell and the Host Response
- Co-stimulation / Immunoregulation
- Dendritic Cells
- Endpoints, Response Criteria for Clinical Trial Design
- Enhancing Cancer Vaccines
- Infectious Agent Vectors
- Innate Immunity to Tumors
- New Agents
- TH-17, Cytokines and T Cell Subsets
- Trafficking and in vivo Imaging
- Tumor Escape / Tumor Microenvironment
- Tumor Targeting Monoclonal Antibodies

Associated Programs

- Global Regulatory Summit - *October 29*
- Primer on Tumor Immunology and Biological Therapy of Cancer - *October 30*
- Workshop on Cancer and Inflammation: Promise for Biological Therapy - *October 30*
- Annual Meeting - *October 31 - November 2*

International Society for
iSBTc
Biological Therapy of Cancer

www.isbtc.org

**“More answers lead to more questions—
repeat as necessary.”**

Timothy Bushnell, PhD Co-Scientific and Technical Director of Flow Cytometry

Read about what inspires
Dr. Bushnell's research at:
bdbiosciences.com/bushnell



Helping all people
live healthy lives

BD supports researchers at every step forward—offering a comprehensive flow cytometry portfolio including instrumentation, reagents, and collaborative services.